



## LOCATOR PAN INSTALLATION

(GX-271/GX-274 Liquid Handlers  
and GX-271/GX-274 ASPEC®)

The locator pan is installed on the locator plate of the instrument. The locator pan holds the tray inserts.

To install the locator pan:

1. Orient the two posts on the bottom of the locator pan toward the back of the instrument.
2. Place the locator pan on the locator plate. The front and right side of the locator pan should be flush with the front and right side of the locator plate and the two posts should be inserted in the locator plate.
3. Locate the two screws included with the locator pan and place them in the holes at the front of the instrument. Using a Phillips screwdriver, tighten the screws.

