Addendum

---

## 506C System Interface User's Guide

Part Number LT3635

This document contains additional information to that provided in the *506C System Interface User's Guide*.

- **Safety** and **Sécurité** sections are added.
- Disregard the **Install the Security Key** section and all other references to the security key.

## Safety

The instrument is intended to be used in a laboratory by trained technical personnel. For safe and correct use, both operating and service personnel should follow the instructions contained in the user's guide when installing, cleaning, and maintaining the instrument.

The following safety precautions must be observed during all phases of operation, service, and repair of the instrument. Failure to comply with these precautions or with specific warnings elsewhere in the user's guide violates safety standards of design, manufacture, and intended use of the instrument. Gilson assumes no liability for the customer's failure to comply with these requirements.

The instrument has been certified to safety standards required in Canada, Europe, and the United States. Refer to the rear panel label on the instrument or the Declaration of Conformity document for the current standards to which the instrument has been found compliant.

The following electronic and hazard symbols may appear on or inside the instrument.

<b>Symbol</b>	<b>Explanation</b>	<b>Symbol</b>	<b>Explanation</b>
~	Alternating current Courant alternatif Wechselstrom		Caution Attention Vorsicht
	Protective conductor terminal Borne de terre de protection Schutzleiteranschluss		Caution, risk of electric shock Attention, risque de choc électrique Vorsicht, Elektroschockgefahr
	Electrical power ON Sous tension Netzschalter ein	○	Electrical power OFF Hors tension Netzschalter aus

## Voltage

Access to the rear panel is necessary. The instrument must be detached from all voltage sources before service, repair, or exchange of parts. For normal operation, the instrument is to be grounded through the AC line cord and power supply provided. Failure to do so can result in a potential shock hazard that could result in serious personal injury.

The instrument must only be operated with the voltage specified on the rear panel label of the instrument and with the grounded AC line cord and power supply provided.

## Replacement Parts

Be sure to use only replacement parts mentioned in the user's guide. Do not repair the instrument or change parts not listed in the user's guide. If it is necessary to replace parts not listed, please contact your local Gilson representative.

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